


ScufaSoft Downloadable Fix

If you receive the following error when trying to start ScufaSoft Software,

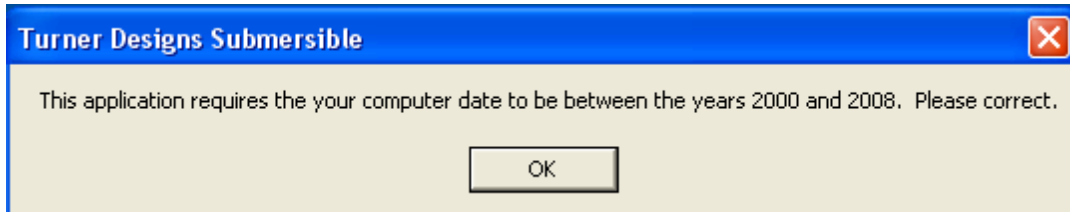

then follow the instructions below, which will help you fix this error.

To download and fix:

- 1) Simply download the ([ScufaSoft.exe.zip](#)) folder to your computer
- 2) Right click on the folder and select 'extract all' to extract the file from the folder
Note: You may use other extracting software if you choose
- 3) Copy the extracted ScufaSoft.exe file to C:\Program Files\Turner Designs\ScufaSoft, which will over-write the existing .exe file
- 4) Then use ScufaSoft as you normally would

If you run into any trouble after the fix has been implemented, please contact Turner Designs' Technical Support at 408-212-4063.